Medix Biochemica

Product Manual Cat. No: #1101

Taq DNA Polymerase Hot-Start

Description

Taq DNA Polymerase Hot-Start is supplied together with the 10x Taq reaction buffer. The reaction buffer has been specifically designed for optimal PCR performance and polymerase activity. Taq DNA Polymerase Hot-Start can also be used for real-time cycling, when adding a suitable real-time dye or a fluorescent probe.

Applications include standard PCR, real-time-PCR (addition of suitable dye or probe required), primer extension reactions, TA cloning, 3'A-tailing of blunt ends, and screening / high-throughput PCRs.

Kit components

Component	S pack*	L pack*
Taq DNA Polymerase Hot-Start	1 x 80 µL	2 x 400 µL
10x Taq reaction buffer	2 x 1.25 mL	13 x 1.25 mL

*Other pack sizes, bulk orders and customization are available upon request.

Storage and shipment

Transport with cool packs. The reagents should be stored at -20°C upon arrival. The reagents are stable until the expiration date if stored correctly.

Reaction master mix set-up

The recommended master mix set-up for a 50 μL reaction volume is shown in the table below.

Reagent	Volume (µL)	Final concentration
Taq DNA Polymerase Hot- Start (5 U/µL)	0.25	1.25 U/rxn
10x Taq reaction buffer	5	1x
∞Forward primer (10 μM)	1	0.2 μM (0.05–1 μM)
∞Reverse primer (10 μM)	1	0.2 μM (0.05–1 μM)
dNTPs (2 nM)	5	200 µM
Template / Sample extract	x	<1000 ng* DNA
Nuclease-free water	Up to 50 µL final volume	

Keep all components on ice.

Spin down and mix all solutions carefully before use.

 ∞ Primers should ideally have a GC content of 40–60% typically.

*Suggested template concentration should be about 1 ng - 1000 ng (genomic DNA) or 1 pg - 1 ng (plasmid/viral DNA) per reaction.

Instrument and program set-up

Cycles	Steps	Temperature	Time
1	Initial denaturation	95°C	2 min
25–40	Denaturation	95°C	15 sec
	Annealing*	54–72°C	30 sec
	Extension	72°C	1 min /1000 bp

*Typically, the annealing temperature is about 3–5°C below the calculated melting temperature of the primers used.



myPOLS Biotec GmbH Byk-Gulden-Strasse 2 78467 Konstanz, Germany VAT no. DE294545185

www.medixbiochemica.com www.mypols.de mdx@medixbiochemica.com

Legal disclaimer

myPOLS Biotec part of Medix Biochemica group products meet their specifications if transported, stored, and used according to the instructions. Medix Biochemica's products may not be used or reproduced without Medix Biochemica's written permission.

17/10/2023 Version 3

Medix Biochemica

Product Manual Cat. No: #1101

Technical information and support

For technical enquiries or assay development support, please contact us via e-mail at: mdx@medixbiochemica.com.

Additional information and technical resources are available on our website at: info.medixbiochemica.com/resources.



myPOLS Biotec GmbH Byk-Gulden-Strasse 2 78467 Konstanz, Germany VAT no. DE294545185 www.medixbiochemica.com www.mypols.de mdx@medixbiochemica.com

Legal disclaimer

myPOLS Biotec part of Medix Biochemica group products meet their specifications if transported, stored, and used according to the instructions. Medix Biochemica's products may not be used or reproduced without Medix Biochemica's written permission.